

# Serum-Free NeoSupp

# #Cat: NB-11-0081 Size: 8ml

### **Product Description:**

NeoSupp Supplement contains bovine growth factor. The 8mL vial activates 500ml Complete Serum-Free Medium, NB-11-0061. Serum-Free NeoSupp is the broad-spectrum mitogen growth supplement designed for use with Serum-Free Medium. Serum-Free NeoSupp contains CSC bovine growth factor and has been designed to supplement growth in the serum-free media environment. NeoSupp is included in various Complete Serum-Free Medium, NB-11-0061 and is also available individually as a stand-alone supplement to improve performance of your serum-free medium.

CSC complete media and reagents are sterile, made with WFI and all components are cGMP and ISO Compliant.

# **Components:**

This 8mL vial supplements 500mL Serum Free Medium. Serum-Free NeoSupp is included in Complete Serum-Free Medium, NB-11-0061.

#### Storage:

Can be frozen at -20°C. Once opened, shelf life is 30 days at +2-8°C. If entire unit will not be used within 30 days, aliquot and freeze.

#### Product Use:

NB-11-0081 is for laboratory research only. It is not approved for human or animal use, or for application in in vitro diagnostic procedures.

# Shipping:

All medium and reagents shipped at ambient temperature.

#### Instructions for Use:

Make sure all work surfaces are disinfected. Spray the outside of your gloves prior to work under the hood.

• Spray the outside of all tools/instruments (bottles, tubes, racks, etc.) with ethanol solution before bringing them under the hood. Use 8mL Serum-Free NeoSupp to supplement 500mL Serum Free cell culture medium, or our Complete Serum-Free Medium, NB-11-0061.

• Open caps on Serum Free medium and NeoSupp, place caps on gauze pads saturated in alcohol.

• Pour or pipette 8mL Serum-Free NeoSupp into 500mL Complete Serum-Free Medium, NB-11-0061.